MarvelXACT™ PEEKsil™

Product Data Sheet



Fused Silica Tubing



Ultra-high



Torque Limiting Technology



Learn More at



100%




Our bioinert MarvelXACT PEEKsil connection systems utilize patented MarvelXACT fittings and PEEKsil tubing for use in the most demanding UHPLC applications. The MarvelXACT fitting features torque-limiting technology that is

finger tight and provides a haptic "click" feedback when it reaches the optimum torque, assuring a perfect zero dead volume connection every time. The robust PEEKsil tubing has smooth silica glass wetted surfaces and is chemically compatible to organic solvents. Each connection system is individually cut, polished, and manufactured to provide industry leading technology that produces superior chromatographic results.

19,000 psi


REUSABLE 100+ USES

FLAT-FACED SEALING CONNECTIONS ZERO DEAD VOLUME

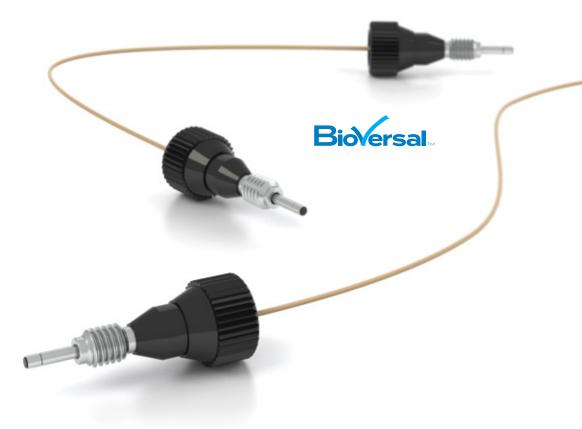





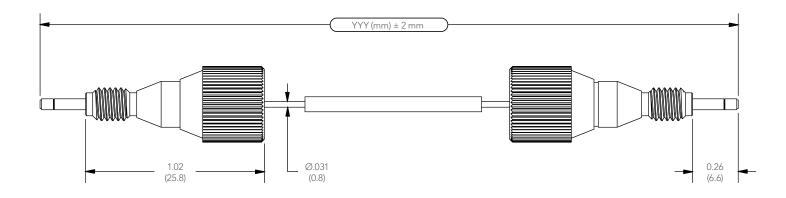

Technical Specifications				
Length:	150 mm	350 mm	600mm	
Bio-inert MarvelXACT PEEKsil Connection Systems				
25 μm ID	UPFF-7025150	UPFF-7025350	UPFF-7025600	
50 μm ID	UPFF-7050150	UPFF-7050350	UPFF-7050600	
75 μm ID	UPFF-7075150	UPFF-7075350	UPFF-7075600	
100 µm ID	UPFF-7100150	UPFF-7100350	UPFF-7100600	

Product Specifications		
Pressure Capability	19,000 psi (~1310 bar, 131 MPa) for Routine Use	
Installation Method	Finger-Tighten Until the First "Click" Feedback is Received	
Tubing Type	1/32" OD PEEK with 1/16" Rigid Tube Ends	
Fitting Type	10-32 Threaded, PEEK Fittings with 316 Stainless Steel Threads	
Wetted Material	PEEKsil	
Max Temp	120°C (Higher Operating Temperatures May Reduce Performance)	
Minimum Tubing Bend Radius	1.0" (~25 mm)	
Max Receiving Port	0.020" (~0.4 mm)	

Note: The above performance specifications apply to use with appropriately designed receiving ports under optimal conditions, using water for the testing process. If different conditions are used, the expected pressure threshold will be different.

IMPORTANT NOTES

MarvelXACT tubing includes a sleeve to assist in product identification, containing details for ID, length, and part number.



REGULATORY COMPLIANCE

As of the date of the publication, MarvelXACT is compliant with current RoHS regulations.

INSTALLATION INSTRUCTIONS

Finger-tighten until the first "click" feedback is received.







Global Premium Distributor:

Biotech AB - Europe | Biotech USA LLC | Bionik Inc. – Japan www.biotechfluidics.com info@biotechfluidics.com | +46 300 56 91 80